

Laboratory Washer

AW62


Model	AW62
Washing Time (Each Process)	0 to 30min. Settable
Washing Water Temp.	45 to 80deg.C

Features :


- All Process (from Washing to Rinsing) are Full Automatic and Each Process are displayed on Indicator
- Washing Process and Time can be set depending on the Container's Shape and Dirty Level
- Washing Water Temp affects washing Result. For this reason, this machine is loaded with Heater to heat up the water.
- Compact Desk Top Type with Turn Table and Powerful Pressurized Water Jet from Two Directions, Top and Bottom.



Dimensions (mm)

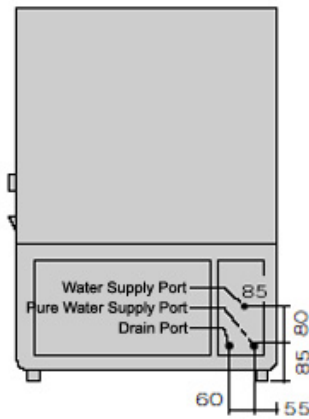

Sequence/Time Chart

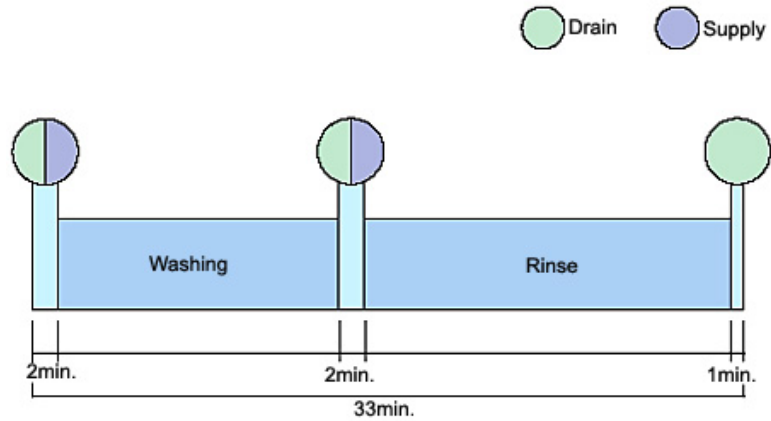

Specifications

Model	AW62
■ Performance	
Washing Type	Tow Direction, Top and Bottom, Pressurized Water Jet Type, Fixed Jet Nozzle (Exchangeable to Jet Rack) Rack Spinning Type
Washing Cycle	Wash (0 to 30min.) Settable Rinse (0 to 30min.) Settable Final Rinse (Option) (Connect with Pure Water Supplier) Washing by Pure Water 20L The above process can be selected by the way of operator's own usage.
Supplying Water Temp.	Room Temp. to 60oC
Washing Temp.	45 to 80deg.C
Hot Water Supply System	Heater Heat-up (Built-in: 6kW)
Supplying Water Pressure	0.1 to 0.3Mpa
Container Stand	Spinning Table (Standard), (Racks are Option)
Water Supplying System	Electromagnetic Valve Open/Close, Water Flow amount is fixed by Liquid Surface Control Switch
Water Drain Type	Natural Drainage by Level Gap
■ Components	
Exterior	Steel Plate, Melamine Resin Paint
Interior	Stainless Steel
External Dimensions	W600 x D620 x H940mm
Inner Dimensions	W560 x D600 x H570mm
Spinning Table	φ550mm (Max. Load 25kg)
Pump	Three-phase, AC200V, 250W
Door	Up/Down, Open/Close Type (Stop at Desired Position)
Weight	Approx. 90kg
Power	Three-phase, AC200V, 50/60Hz, 30A
■ Accessory	
Supply Hose (with Coupler)	2m, 1pc.
Drain Hose	I.D. 25.4mm, 2m, 1pc.
Vinyl Cover	
Amorphous Phosphorus Detergent	1kg
Main Jet Nozzle Washing Brush	1pc.
Water Supply Port Unit	

Optional Item (for AW62)

Pure Water Supplier	With 20L Pure Water Tank
Jet Rack	42pcs. Settable in case of 100ml Volumetric Flask
Beaker Rack	Approx. 85pcs. settable in case of 50ml Beaker
Test Tube Rack	Approx. 600pcs. settable in case of Φ 16.5mm Test Tube
Flask Rack	Approx. 68pcs. settable in case of 60ml Flask
Amorphous Phosphorus Detergent	8kg



Pure Water Supplier



AW62+Pure Water Supplier



Jet Rack



Beaker Rack



Test Tube Rack



Flask Rack



Amorphous Phosphorus Detergent